

Wavelength Meter

Next Generation

WR8-2



HighFinesse
The Standard of Accuracy



Ångstrom

Available Measurement Range

WR8-2⁹⁾ Standard 330 – 1180 nm

Absolute (and Other) Accuracies¹⁾

330 – 375 nm 10 MHz^{2) 3)}

375 – 1180 nm 2 MHz^{2) 3)}

Quick coupling accuracy (with 50 µm multi mode fiber) 100 MHz

Wavelength deviation sensitivity/Measurement resolution⁴⁾ 0.1 MHz⁵⁾

Measurement Speed⁷⁾

1000 Hz

Required Minimum Input Energy and Power⁸⁾

Standard (VIS) 0.08 – 60 µJ or µW

FSR of the Fizeau Interferometers (Fine/Wide Mode)

2 GHz/20 GHz⁶⁾

1) According to 3σ criterion, but never better than 20 % of the laser linewidth.

2) ± 2 nm around calibration wavelength. Outside of this range, the accuracy as WR8-10.

3) ± 200 nm around calibration wavelength. Outside of this range, the accuracy as WS7-30.

4) Standard deviation. WR8-2 requires photonic crystal (endlessly singlemode) fibers to reach this resolution.


5) 100 kHz for special ranges on request.

6) Each instrument in each mode can measure lasers with a linewidth up to 30 % of the corresponding FSR.

7) Depending on PC hardware and settings.

8) The CW power interpretation in [µW] compares to an exposure of 1s (generally the energy needs to be divided by the exposure time to obtain the required power).

9) Next Generation (NG).

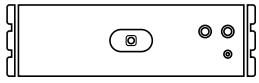

Wavelength Meter
Next Generation

WR8-2



HighFinesse
The Standard of Accuracy



Ångstrom

Calibration

Recommended calibration period ≤ 2 minutes


SLR-780 Rack or any well known laser source $\Delta\nu < 1$ MHz

Warm-up Time

> 30 min. warm-up, or until ambient equilibrium

Dimensions L \times W \times H¹⁰⁾

541 \times 483 \times 133 mm

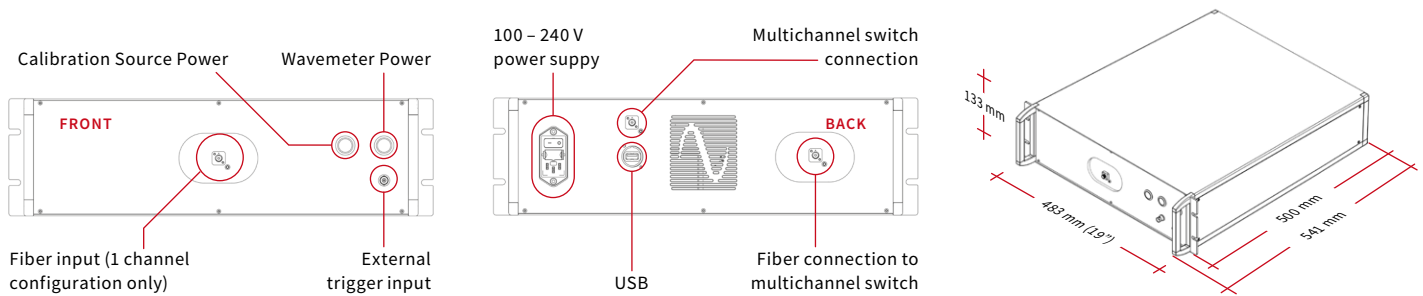

Weight

16 kg

Interface


Control High-speed USB 2.0 connection

External Trigger BNC

Power Supply

100 – 240 V ~ 50/60 Hz 80 W

10) Dimensions with handles.

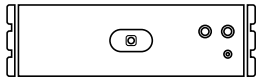

Wavelength Meter
Next Generation

WR8-2



HighFinesse
The Standard of Accuracy



Ångström

Included Options

External Trigger (TTL)

The TTL option guarantees exact synchronization of the wavelength measurement with other measurement tasks in your lab. It provides low-noise signals without parasitic parts when measuring pulsed signals with low duty cycles. The wavelength measurement can be triggered by sending a standard 5 V TTL signal to the input of the wavemeter (for example generated by a delay generator). Necessary cables to connect to a BNC output are included.

Please note, if the option MC is ordered together with the TTL option, the TTL mode can only be used if the switch is set fixed to one input channel.

Photonic Crystal Switch Rack

In order to measure the frequencies of more than just one laser at a time, an opto-mechanical switch is used. The combination of our high-speed wavelength meters with one of the quickest fiber switches available allows up to eight channels to be measured almost simultaneously. Exposure time and other parameters can be defined independently for each light source.

The WR8-2 series features the use of an endlessly singlemode switch based on the photonic crystal switch technology in an external 19", 1 HU rack. This allows to measure any laser wavelength on all switch input channels within all measurement ranges of the WR8-2.

The TTL mode can only be used if the switch is set fixed to one input channel.

Options

Laser Control (PID)


With the PID option it is possible to stabilize the frequency of a laser connected to the wavelength meter using a software based proportional-integral-derivative controller (PID controller). Unlike analog PID electronics, the PID option provides software based signal processing, allowing the laser to be stabilized to a specific user defined frequency or regulated with an arbitrary pattern.

This makes it extremely useful in experiments where the laser frequency has to be actively regulated or varied to fit changing experimental conditions, such as laser cooling, atomic detection, trapping and spectroscopy.

Combined with the MC option the wavelength meter can be used to stabilize multiple lasers simultaneously. The regulation speed, quality and absolute accuracy match the measurement speed, relative accuracy and absolute accuracy of the wavelength meter respectively. The measurement speed is not affected by the regulation.

Multichannel Laser Control (MCLC)

The MCLC option combines the capabilities of the photonic crystal switch and PID option in one instrument. It can be ordered with 4 or 8 channels. This solution for 4 or 8 channel PID is compact and circumvents the need for a computer with a free PCIe slot such that it is possible to monitor and control up to 8 lasers using the HighFinesse wavemeter connected to a laptop as the 4 or 8 analogue outputs with maximum ± 10 V are located at the MCLC instrument. At the same time the MCLC guarantees for the same excellent feedback rate and precision the HighFinesse PID option is known for. Adapters to connect from the MCLC to the laser from Lemo 00 to BNC are included in the scope of delivery. The MCLC is available as a tabletop system only.

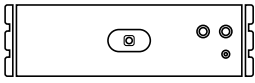

Wavelength Meter
Next Generation

WR8-2



HighFinesse
The Standard of Accuracy



Ångström

Options

External Calibration Sources for Rack Systems

Our rack calibration sources are highly accurate, reliable, plug and play and ready to be mounted in a 19" Rack system. This guarantees the accuracy and stability of measurements with our wavelength meters. We offer a variety of frequency stabilized, narrow linewidth, laser sources with up to ± 0.5 MHz frequency stability for different applications.

For further information see our product description here: <https://www.highfinesse.de/cal>

Typical Applications

The WR8-2 is the highest-end solution for wavelength monitoring and control with an absolute accuracy of 2 MHz and a wavelength deviation sensitivity of 0.1 MHz. It comes with the photonic crystal fiber technology enabling multichannel operation in the spectral range of the wavelength meter.

The 2 MHz accuracy presumes, that the instrument is supported by one of our calibration sources based on atomic spectroscopy (or an laser source provided with comparable accuracy).

Further Information

For further technical information, application examples, diagrams and for customization of the WR8-2 please contact:

HighFinesse Team

service@highfinesse.de



HighFinesse GmbH
Wöhrdstraße 4
72072 Tübingen, Germany



T +49 (0) 7071 - 53918 0
F +49 (0) 7071 - 53918 99
M info@highfinesse.com



Additional information
and distributors:
www.highfinesse.com

